



Operating the Wig-L-Bug with 1" long vials (plastic or steel)

The #3113 adapter is used with 1" long vials (plastic or steel). The black anodized adapter is mounted on the arms as follows: The long black side across the top of the two arms and the two short black sides on the outside of the arms facing down. The biggest screw outside the front arm and the small screw and nut for the rear arm. Loosen screw and nut completely. Place the short screw in hole of the rear arm clear thru the back slotted side of the adapter (adapter flange on the outside of arm). Put on nut and tighten. Adapter is now loose at the front. Place the 1" vial between the arms and place the adapter, long side over top of arms - front short black flange outside the front arm. Holes on top of adapter should slip over arms slightly. Then tighten the front (10/32) socket screw to hold vial steady.